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Home

Search & Browse for siRNAs & Assays

Search Database to Order siRNAs & Assays
Search by gene name, gene symbol, gene transcript ID, or product ID. [Search Database >](#)

Browse Pathways to Order siRNAs & Assays
Browse to cellular pathway diagram, then click protein to see associated products. Also download full product list for pathway. [Browse Pathways >](#)

Order Custom siRNA

Order Custom Designed siRNA
When your needs extend beyond our vast collection of pre-designed siRNAs ... [Order Custom >](#)

Order Custom Synthesized siRNA
siRNAs synthesized using sequences you provide. [Order Custom >](#)

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Next

Choose Application (Product to Search)

- For use in cells (Silencer® Select siRNAs)
- For use in animals (Ambion® In Vivo siRNAs)

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Search Database to Order siRNAs & Assays

Recent Searches:

Field to be Searched	Criteria	Species	Search Term	
NCBI Transcript	Contains	All Species	NM_000586	<input type="submit" value="Submit"/>

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Search Database to Order siRNAs & Assays

Recent Searches: [NM_000586](#)

Field to be Searched	Criteria	Species	Search Term	
<input type="text" value="NCBI Transcript"/>	<input type="text" value="Contains"/>	<input type="text" value="All Species"/>	<input type="text" value="NM_000586"/>	<input type="button" value="Submit"/>

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Search Database to Order siRNAs & Assays

Recent Searches: [NM_000586](#)

Field to be Searched	Criteria	Species	Search Term	
NCBI Transcript	Contains	All Species	NM_000586	Submit

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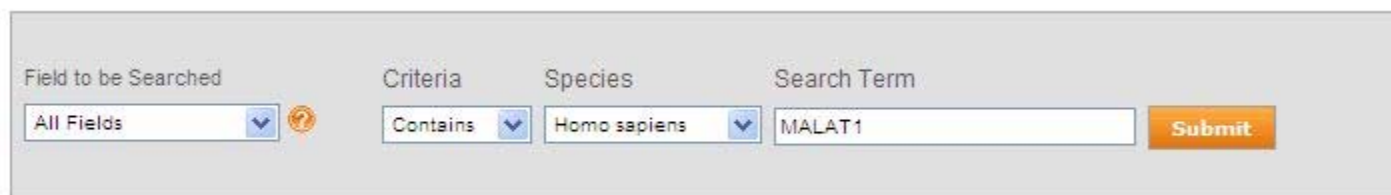
siRNAの検索 Noncoding Transcriptの検索

Field to be Searched: All Fields
Criteria: Contains
Species: Homo sapiens
Search Term: MALAT1
Submit

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The screenshot shows a web interface for searching siRNAs. At the top, there is a navigation bar with buttons for Home, Search, Select siRNAs/Assays, Select Other Products, and Review & O. Below this, there is a section titled "Choose Application (Product to Search)" with two radio buttons: "For use in cells (Silencer[®] Select siRNAs)" (selected) and "For use in animals (Ambion[®] In Vivo siRNAs)". A "More Information" link is also present. The next section is "Search Database to Order siRNAs & Assays", which includes a "Recent Searches" list with "NM_000586". Below this is a search form with the following fields: "Field to be Searched" (set to "NCBI Transcript"), "Criteria" (set to "Contains"), "Species" (set to "All Species"), and "Search Term" (set to "NM_000586"). A "Submit" button is located to the right of the search term field. Below the search form, the "Search Results" section displays "1 result(s) for NCBI Transcript contains 'NM_000586'". The results are shown in a table with the following columns: "NCBI Transcript", "Gene ID / RNAdb ID", "Gene Symbol / RNAdb Gene Symbol", "Gene Name / Description", "Species", and a "Select" button. The "Select" button for the first result is highlighted with a red box.

NCBI Transcript	Gene ID / RNAdb ID	Gene Symbol / RNAdb Gene Symbol	Gene Name / Description	Species	
NM_000586.3	3558 / -	IL2 / -	interleukin 2 / -	Homo sapiens	Select

siRNAの検索 V

Select siRNAs & Assays

Gene ID / RNAdb ID	Gene Symbol / RNAdb Gene Symbol	Gene Name / Description	Species	Product Count
3558 / -	IL2 / -	interleukin 2 / -	Homo sapiens	29
Gene Alias:	IL-2 TCGF lymphokine			
Transcripts:	NM_000586 Additional Information			
Pathways:	View 70 pathways			

[Alignment Map](#)

Unable to find the siRNAs you need? Order [Custom Designed siRNAs](#).

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Home Search **Select siRNAs/Assays** Select Other Products Review & Order

← Previous Next →

Select siRNAs & Assays

Gene ID / RNAdb ID	Gene Symbol / RNAdb Gene Symbol	Gene Name / Description	Species	Product Count
3558 / -	IL2 / IL2	Interleukin 2 / -	Homo sapiens	29

Gene Alias: IL-2 TCGF lymphokine
Transcripts: [NM_000586 Additional Information](#)
Pathways: [View 70 pathways](#)

Alignment Map
Unable to find the siRNAs you need? Order [Custom Designed siRNAs](#).
Silencer® Select siRNAs For Coding Transcripts [\[More Info\]](#)

siRNA ID	Type	Availability	Size/Purity	Quantity	Add All
s223931	Pre-designed: Coding	Inventoried	5 nmol / STD: 4392420	1	Add
s7283	Pre-designed: Coding	Inventoried	5 nmol / STD: 4392420	1	Add
s223932	Pre-designed: Coding	Inventoried	5 nmol / STD: 4392420	1	Add

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Home Search Select siRNAs/ Assays Select Other Products Review & Order

Please review your Shopping List below and make any desired modifications, then click "Add to Basket" to begin the purchasing process.

If not already logged in, you will be prompted to log in to the Applied Biosystems Store before adding products to basket. If you do not have a Store account, registration is quick and easy. Register with AB Store

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Select All 製品

! You must order at least 4 siRNAs at the 1 nmol scale. Please add additional siRNAs or change the scale to 5 nmol.

1. Silencer® Select Pre-designed siRNA, inventoried, std purity, 1 nmol

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Select Other Products

Review & Order

Please review your Shopping List below and make any desired modifications, then click "Add to Basket" to begin the purchasing process.

If not already logged in, you will be prompted to log in to the Applied Biosystems Store before adding products to basket. If you do not have a Store account, registration is quick and easy. [Register with AB Store](#)

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Store Log In

Use this page to register and/or login to the Applied Biosystems Store. If you have [Us](#).

Store Log In

To login to the Applied Biosystems Store, enter your Username and Password below, and then click **Login**.

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Submit Your Shopping List

Select a shopping basket or create a new one for your assays. Then click the Add to Basket button.

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Organization	Life Technologies Japan Ltd.	PO Box Postal Code	
Department		City	Chuo-ku,
Attention		District	Tokyo
Building/Room		Postal Code	1040032

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The screenshot shows the top navigation area of the Applied Biosystems website. On the left is the Applied Biosystems logo. To the right are links for 'Store Log Out', 'My Account', 'Shopping Basket', and 'Order Inquiry'. A search bar is also present. Below these is a main navigation bar with 'Home', 'Products & Services', 'Support', and 'About Us'. A secondary navigation bar contains 'Quick Order', 'Favorites', 'Workspace', 'Manage Shopping Baskets', 'Promotions/Quotes', and 'Store Help'. At the bottom of the navigation area are numbered steps: 1 Shopping Basket, 2 Shipping & Billing, 3 Summary, and 4 Confirmation.

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Number of Designs: (Maximum 6)

Apply *Silencer*® Select siRNA chemical modifications? Yes No ← siRNAの種類を選択

Silencer® Select siRNA chemical modifications reduce off-target effects; they do not diminish silencing efficiency.

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Custom Designed Silencer® Select siRNAs

siRNA ID	Type	Availability	Purity/Size	Quantity	Add All
ABGJPBZ	Custom	Made to Order	4399889: HPLC 40 nmol	1	Add
ABHSNH7	Custom	Made to Order	4399889: HPLC 40 nmol	1	Add
ABI1LOF	Custom	Made to Order	4399889: HPLC 40 nmol	1	Add

Also available... these Pre-Designed/Validated siRNAs that target your gene/transcript of interest.

Silencer® Select siRNAs [More Info]

siRNA ID	Type	Availability	Purity/Size	Quantity	Add All
s7327	Validated	Inventoried	4390824: STD 5 nmol	1	Add
- Validated by assay Hs00174103_m1					
s223934	Pre-Designed	Inventoried	4392420: STD 5 nmol	1	Add
s7328	Validated	Inventoried	4390824: STD 5 nmol	1	Add
- Validated by assay Hs00174103_m1					

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Shopping List (0 items)
Order Now Clear List

Custom Designed Silencer® Select siRNAs

siRNA ID	Type	Availability	Purity/Size	Quantity	Add All
ABGJPBZ	Custom	Made to Order	4399889: HPLC 40 nmol	1	Add
ABHSNH7	Custom	Made to Order	4399889: HPLC 40 nmol	1	Add
ABI1LOF	Custom	Made to Order	4399889: HPLC 40 nmol	1	Add

Also available... these Pre-Designed/Validated siRNAs that target your gene/transcript of interest.

Silencer® Select siRNAs [More Info]

siRNA ID	Type	Availability	Purity/Size	Quantity	Add All
s7327	Validated	Inventoried	4390824: STD 5 nmol	1	Add
- Validated by assay <i>Hs00174103_m1</i>					
s223934	Pre-Designed	Inventoried	4392420: STD 5 nmol	1	Add
s7328	Validated	Inventoried	4390824: STD 5 nmol	1	Add
- Validated by assay <i>Hs00174103_m1</i>					

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Home > **Input Design** > Confirm Design > Select Other Products > Review & Order

Next →

Order a Custom Designed siRNA

Name Your Target

This name will help identify your siRNA(s) during order processing and will be printed on the label. We recommend using the gene symbol when possible.

Target Name: (8 characters max.)

Provide Target Information

Target Format: ▼

Entrez ID:

Select Options

Number of Designs: (Maximum 6)

Apply *Silencer*[®] Select siRNA chemical modifications? Yes No

Silencer[®] Select siRNA chemical modifications reduce off-target effects; they do not diminish silencing efficiency.

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Home > **Input Design** > Confirm Design > Select Other Products > Review & Order

Next → Shopping List (0 items)

Order a Custom Designed siRNA

Advanced Specificity Options

I would like an additional off-target check performed against this species: ←

e.g. You would choose "mouse" if you are targeting the transcripts of a human gene and, in addition to not targeting any other human transcripts, you would like the siRNA to not target any mouse transcripts (except for any specified to be targeted below).

I would like this siRNA to also target not target this

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Home > Input Design > **Confirm Design** > Select Other Products > Review & Order

Order a Custom Designed siRNA

Target Name: Testsi
Select Modifications: Yes
Target Format: Entrez Gene ID
Target Entrez Gene ID: 3576

[Edit](#)

Custom Designed *Silencer*[®] Select siRNAs

siRNA ID	Type	Availability	Purity/Size	Quantity	Add All
ABGJPBZ	Custom	Made to Order	4399885: STD 5 nmol	1	Add
ABHSNH7	Custom	Made to Order	4399885: STD 5 nmol	1	Add
ABI1LOF	Custom	Made to Order	4399885: STD 5 nmol	1	Add

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GeneAssist™ siRNA Workflow Builder

Home

Search & Browse for siRNAs & Assays

Search Database to Order siRNAs & Assays

Search by gene name, gene symbol, gene transcript ID, or product ID.

Search Database >

Browse Pathways to Order siRNAs & Assays

Browse to cellular pathway diagram, then click protein to see associated products. Also download full product list for pathway.

Browse Pathways >

Products

- *Silencer*® Select Pre-designed & Validated siRNAs
- *Silencer*® Pre-designed & Validated siRNAs
- TaqMan® Gene Expression Assays

Order Custom siRNA

Order Custom Designed siRNA

When your needs extend beyond our vast collection of pre-designed siRNAs ...

Order Custom >

Order Custom Synthesized siRNA

siRNAs synthesized using sequences you provide.

Order Custom >

Products

- Custom Designed *Silencer*® Select siRNA
- Custom Designed *Silencer*® siRNA
- Custom Select siRNA
- Custom siRNA

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GeneAssist™ siRNA Workflow Builder

Home > **Input Sequence** > Confirm Sequence > Select Other Products > Review & Order

Next →

Order a Custom Synthesized siRNA

Select Quantity and Format

Number of siRNAs to be synthesized: (Maximum 50)

Sequence Format: siRNA Sequence ← Target (mRNA) Sequence

Select Options

Apply *Silencer*® Select chemical modifications? Yes No

Silencer® Select chemical modifications reduce off-target effects; they do not diminish silencing efficiency.

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GeneAssist™ siRNA Workflow Builder

Home > **Input Sequence** > Confirm Sequence > Select Other Products > Review & Order

Order a Custom Synthesized siRNA Next →

Enter Sequence Information [Silencer³ Select siRNA Modification On]

siRNA Name: Purity/Size: 4390828: STD 20 nmol

Sense Sequence: (Passenger Strand) 5' 3' (21 nt) Must be 21 nt long.

Antisense Sequence: (Guide Strand) 5' 3' (21 nt)

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GeneAssist™ siRNA Workflow Builder

Home > Input Sequence > **Confirm Sequence** > Select Other Products > Review & Order

Next ➔

Order a Custom Synthesized siRNA

Sequence Format: siRNA Sequence
Select Modifications: Yes

Custom Synthesized *Silencer*® siRNAs

siRNA ID	Type	Availability	Part #, Purity, Size	Quantity	Add All
ADBJWLC	Custom	Made to Order	4390828: STD 20 nmol ▼	1	Add
Edit	siRNA Name: target01 Sense: 5' GACUGAUCGAGUCGAGUCUITT 3' Antisense: 3' GACUGACUAGCUCAGCUCAGA 5' Special Instructions:				

Shopping List (0 items)

Order Now Clear List

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GeneAssist™ siRNA Workflow Builder

Home **Input Sequence** Confirm Sequence Select Other Products Review & Order

Next →

Order a Custom Synthesized siRNA

Select Quantity and Format

Number of siRNAs to be synthesized: (Maximum 50)

Sequence Format: siRNA Sequence Target (mRNA) Sequence

Select Options

Apply *Silencer*® Select chemical modifications? Yes No

Silencer® Select chemical modifications reduce off-target effects; they do not diminish silencing efficiency.

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Next →

Order a Custom Synthesized siRNA

Sequence Format: Target (mRNA) Sequence
Select Modifications: Yes

Custom Synthesized *Silencer*® siRNAs

siRNA ID	Type	Availability	Part #, Purity, Size	Quantity	Add All
ADCSURK	Custom	Made to Order	4390828: STD 20 nmol	1	Add

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The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners.

TaqMan is a registered trademark of Roche Molecular Systems, Inc

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©2010 Life Technologies Japan Ltd. All right reserved.